

PREMIER Biosoft

Accelerating Research in Life Sciences





How to import AB SCIEX's .wiff file?



SimGlycan[®]

Step 1: Install and activate SimGlycan

- Install the program by double clicking the appropriate installer.
- After installation, activate the program using your registration number and email address.

Biosoft

• Once registration is complete, close the program.

Step 2: Check if .NET framework version (v 4.0.30319) is installed

18 I 📑 18 🖛 I			v4.0.303	19		×
File Home Share V	ïew				~ ~	?
🛞 🍥 👻 🕇 퉬 > This PC	▹ OS (C:)	t.NET → Framework → v4.0.30319	•		✓ C Search v4.0.30319	ρ
🔆 Favorites	Name	Date modified	Туре	Size		^
🔜 Desktop	퉬 1033	3/18/2014 2:55 PM	File folder			
🗼 Downloads	퉬 ASP.NETWebAdminFiles	8/22/2013 9:06 PM	File folder			
📳 Recent places	퉬 Config	8/22/2013 9:06 PM	File folder	G	o to:	
	퉬 en-US	3/18/2014 2:55 PM	File folder			
	퉬 MSBuild	8/22/2013 9:06 PM	File folder	'C	:\Windows\Microsoft.NEI\Framework	
	遇 MUI	8/22/2013 9:06 PM	File folder		A 0 30319 ⁽	
	퉬 Nativelmages	3/18/2014 3:36 PM	File folder		4.0.00010	
	퉬 SQL	8/22/2013 9:06 PM	File folder			
	퉬 WPF	3/18/2014 3:36 PM	File folder	lf f	this folder doesn't exist then download	
🖏 Homegroup	🚳 Accessibility.dll	8/10/2013 6:21 AM	Application extens	36 KE		
	AddInProcess.exe	8/10/2013 6:21 AM	Application	30	nd install the .NET Framework v	
💻 This PC	AddInProcess.exe.config	6/18/2013 5:57 PM	CONFIG File	1 KF	n 20240	
Ҏ dell (badri)	AddInProcess32.exe	8/10/2013 6:21 AM	Application	42 KE	0.30313	
Lesktop	AddInProcess32.exe.config	6/18/2013 5:58 PM	CONFIG File	1 KE	В	
Documents	AddlnUtil.exe	8/10/2013 6:21 AM	Application	42 KE	В	
Downloads	AddInUtil.exe.config	6/18/2013 5:58 PM	CONFIG File	1 KE	В	
Music	🚳 AdoNetDiag.dll	8/10/2013 6:21 AM	Application extens	167 KE	В	
Pictures	adonetdiag.mof	6/18/2013 5:58 PM	MOF File	8 KE	В	
Videos	adonetdiag.mof.uninstall	6/18/2013 5:58 PM	UNINSTALL File	2 KE	В	
L OS (C:)	🚳 alink.dll	8/10/2013 6:21 AM	Application extens	115 KE	В	
New Volume (D:)	AppLaunch.exe	8/10/2013 6:22 AM	Application	97 KE	В	
	applaunch.exe.config	8/22/2013 9:04 PM	CONFIG File	1 KE	В	
📬 Network	Aspnet.config	8/22/2013 9:04 PM	CONFIG File	1 KE	В	
т	aspnet_compiler.exe	8/10/2013 6:22 AM	Application	55 KE	В	
	aspnet_filter.dll	8/10/2013 6:22 AM	Application extens	34 KE	В	
	🚳 aspnet_isapi.dll	8/10/2013 6:22 AM	Application extens	25 KE	В	
	🚳 Aspnet_perf.dll	8/10/2013 6:22 AM	Application extens	41 KE	В	
	aspnet_perf.h	6/18/2013 5:58 PM	H File	8 KE	В	~
349 items						

PREMIER Biosoft

Step 3: Download the required Microsoft .NET framework version

Install it using "dotNetFx40_Full_x86_x64.exe" by downloading "Microsoft .NET Framework 4 (Standalone Installer)" directly from Microsoft website at:

http://www.microsoft.com/en-in/download/details.aspx?id=17718

Biosoft

SimGlycan[®]

Step 4: Launch Command Prompt with Administrator privilege





Launch Command Prompt as Administrator (for Win 7)

ER Biosoft

Type **cmd** in the search field. Upon suggesting **cmd.exe**, right click on it and press **Run as administrator**.

Programs (1)		Open
cmd eve	۲	Run as administrator
Documents (11)		Open with MSConvertGUI
		Open with SeeMS
libtabix.a		7-Zip
📄 libbam.a		Edit with Notepad++
knetfile.h	6	Scan with Microsoft Security Essentials
knetfile.h		Add to archive
changelog.txt		Add to "cmd.rar"
🔄 libtabix.a		Compress and email
libbam.a		Compress to "cmd.rar" and email
Files (170)		Pin to Taskbar
Cmd.pm		Pin to Start Menu
Cmd.pm		Restore previous versions
mk_cmds.1.gz		Send to
Caffeine1 - PubChem.htm		Cut
g pubchem.widgets.js		Сору
S core156.js		Paste
Crmd		Delete
		Open file location
		Properties

Step 5: Execute the first command

• Copy the following command:

cd "%Appdata%\PBI\SGEE\NativeLibs"

 Go to Command Prompt, right click on it, press Paste and then press Enter.

Microsoft Window	Administrator: Command Prompt – – ×
(c) 2013 Microso C:\Windows\syste	t Corporation. Hil Fights reserved.
	Mark Copy Enter Paste
	Select All Scroll Find
and the state of t	PREMIER Biosoft

Step 6: Execute the second command

- Similarly, run the second command
- C:\Windows\Microsoft.NET\Framework\v4.0.30319\RegAsm.exe /codebase WiffReaderCOMWrapper.dll /tlb

	Administrator: Command Prompt - 🗆 🗙	
	Microsoft Windows [Version 10.0.16299.309] (c) 2017 Microsoft Corporation. All rights reserved.	<u>^</u>
	C:\WINDOWS\system32>cd "%Appdata%\PBI\SGEE\NativeLibs"	
	C:\Users\Admin\AppData\Roaming\PBI\SGEE\Nativelibs> <mark>C:\Windows\Microsoft.NET\Framework\v4.0.30319\RegAsm.e</mark> xe /codebase WiffReaderCOMWrapper.dll /tlb	\bigcirc
		FD Biosoft
10.0		V DIUSOII

Step 7: Verify the execution status

The Execution console output should display 'Types registered successfully.'



ER Biosoft

This message indicates that SimGlycan is ready for importing .wiff files.

Step 8: Import .wiff file in SimGlycan

3	SimGlycan® Enterprise Client 5.2 - C:\Program Files\SimGlycan Enterprise Client 5.2_Beta\DemoProject\Wiff_import.sgp – 🗇 🛛 🛛				
File Edit View Analyze Online Tools Help	p				
🚡 🖾 🖉 🕬 🕬	***** D = > * ‱ * * * * * * * * * * * * * * * * *				
Project Navigator ×					
Wiff import Vinf import Vinf import Vinf import Vinf import Sequence Draw Glycan Glycopeptide LC-MS and MS/MS Data Glycan Quantitation	Open MS/MS Data Files × Look in: AB SCIEX Image: Control of the second s				
	Recent Items Desktop Documents This PC Network				
	File name: Open				
	Files of type: WIFF Files (*.wiff) Cancel				
	Load intensity: C Percent (Relatively) 💿 Absolute Value > 🔽 0 %				
ppens an Ab SCIEX Data Tile		ja			
	PRE 0	MIER Biosoft			

Thank You!

2

PREN

MIER Biosoft